

## KAPA SYBR® FAST One-Step qRT-PCR Kits

*One Evolved Polymerase. One Step. One Tube.*

**KAPA SYBR FAST One-Step qRT-PCR Kits** are designed for optimal performance in stringent qRT-PCR reaction conditions, exhibiting dramatic improvements in sensitivity, specificity, reaction efficiency, and fluorescence. The kit is optimized for rapid one-step, one-tube RNA quantification, reducing experimental variation and contamination with a convenient qRT-PCR protocol.




The kits contain M-MuLV reverse transcriptase, RNase inhibitor, and a novel DNA polymerase evolved via directed evolution.

Kits offer:

- SNP genotyping
- Gene expression analysis
- Microarray validation
- Gene knockdown validation
- RNAi and miRNA research

### Increase signal and reaction efficiency




- Consistently higher amplification efficiencies for accurate qPCR
- Higher fluorescence, earlier Cq values and improved reaction efficiencies

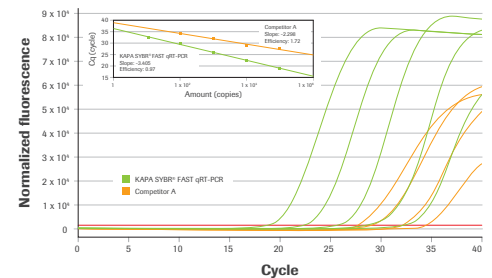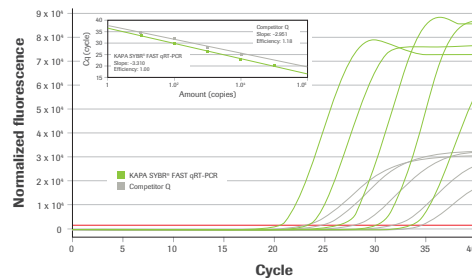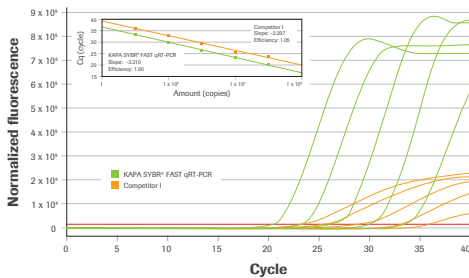

RRM1 gene (94 bp, 45.7% GC) amplified from 10-fold dilution series of RNA (100 ng to 10 pg per 20  $\mu$ L reaction) isolated from human placenta. KAPA SYBR FAST One-Step qRT-PCR Kit demonstrates earlier Cq values, higher fluorescence, and optimal reaction efficiencies.

## Amplifier across a broad range of target lengths


- Robust performance independent of amplicon size

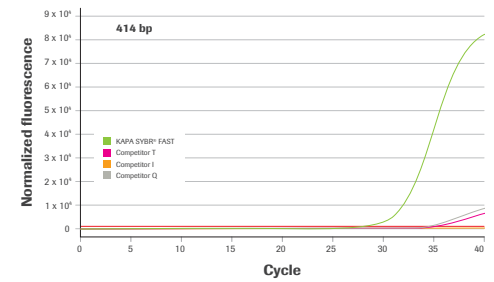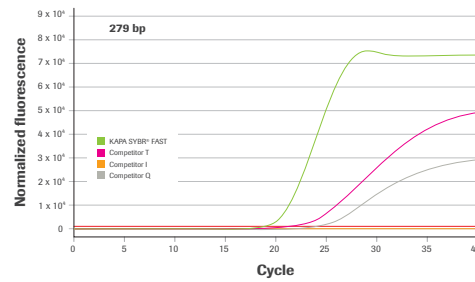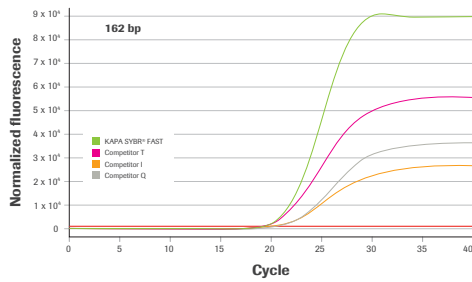

MS2 bacteriophage RNA targets of increasing length were amplified at 0.8 pg per 20  $\mu$ L reaction. KAPA SYBR<sup>®</sup> FAST One-Step qRT-PCR Kit maintained high performance across all amplicon lengths.

## Improve sensitivity and reproducibility

- Accurate interrogation across a wide range of RNA template concentrations

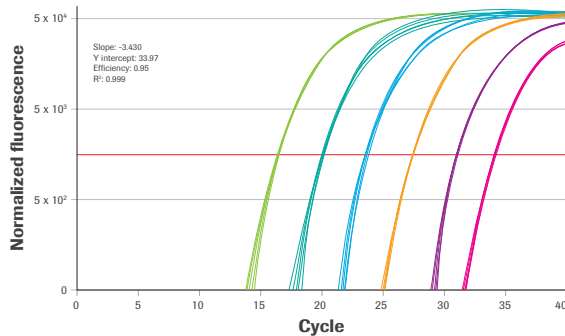

279 bp fragment amplified in replicates of 6 from a 10-fold dilution series of commercial MS2 bacteriophage RNA (80 pg to 0.8 fg per 20  $\mu$ L reaction). Reaction efficiency of 95% calculated for all 6 replicates per data.

## Ordering information

Roche cat. no.	KAPA code	Description	Kit size
07959613001	KK4650	KAPA SYBR <sup>®</sup> FAST One-Step Universal	1 x 1 mL
07959621001	KK4651	KAPA SYBR FAST One-Step Universal	1 x 5 mL
07959630001	KK4652	KAPA SYBR FAST One-Step Universal	2 x 5 mL
07959656001	KK4660	KAPA SYBR FAST One-Step ABI Prism <sup>®</sup>	1 x 1 mL
07959664001	KK4661	KAPA SYBR FAST One-Step ABI Prism	1 x 5 mL
07959672001	KK4662	KAPA SYBR FAST One-Step ABI Prism	2 x 5 mL
07959702001	KK4670	KAPA SYBR FAST One-Step Bio-Rad iCycler <sup>™</sup>	1 x 1 mL
07959729001	KK4671	KAPA SYBR FAST One-Step Bio-Rad iCycler	1 x 5 mL
07959737001	KK4672	KAPA SYBR FAST One-Step Bio-Rad iCycler	2 x 5 mL
07959753001	KK4680	KAPA SYBR FAST One-Step for Roche LightCycler <sup>®</sup> 480	1 x 1 mL
07959761001	KK4681	KAPA SYBR FAST One-Step for Roche LightCycler <sup>®</sup> 480	1 x 5 mL

Published by:

### Roche Sequencing and Life Science

9115 Hague Road  
Indianapolis, IN 46256

[sequencing.roche.com](http://sequencing.roche.com)

[go.roche.com/SybrFastOneStep](http://go.roche.com/SybrFastOneStep)

For Research Use Only. Not for use in diagnostic procedures.

KAPA, LIGHTCYCLER, and SYBR are trademarks of Roche. All other product names and trademarks are the property of their respective owners.

© 2020 Roche Sequencing and Life Science. All rights reserved.

MC-US-06989

SS202001

A520

6/20